

# Release Notes

---

## Internationalization Note

- For release 2.2, this software is **only** recommended and validated while running with US-English default regional settings. Running with other regional settings is not supported.

## Envoy 2.2.3.1

---

In this release, there is an additional install shortcut to run Hydra IIc VHC. To be able to run this customized method process, the software will require a License Dongle to be attached to the controlling computer.

## New Features added:

### Installation of Communication Drivers

---

The Installation of the USB drivers that facilitate communication between the software and hardware has been greatly simplified and incorporated into the software installation of Envoy 2.2.

### Hydra IIc Sleep Button Behavior

---

Cooling the catalyst with Oxygen flowing and assuring the catalyst is sealed from room air when the instrument will be inactive has been automated.

### Significant Digits on Calibration Display and Printout

---

The ability to customize the number of significant digits (or the number of digits after the decimal point) in the Calibration Statistics Table below the Calibration Graph has been added.

Additionally, on the Results and Report tabs, the “Hg [ng]” and the “Weight [g]” columns are now controlled by the user selected menu item “Analysis Report Precision”.

### Calibration Curve Display

---

The X and Y axes and the Plot of the Calibration Graph have been reformatted to align with the calibration equation.